


# S.NEST 2.0 C.NEST 2.0 C.BIRD

Early Stage Mixing at 24/96 Well  
Scale — Built for Clone Selection

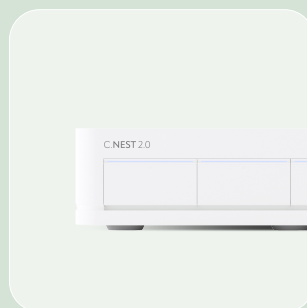


The Leadgene microbioreactor family, including S.NEST<sup>®</sup>, C.NEST<sup>®</sup>, and C.BIRD<sup>™</sup>, forms a unified bioprocessing platform that supports cell line development from single cell expansion through early stage screening.

Built on three core technologies, programmable mixing in 96 and 24 well plates, real time DO and pH monitoring, and four independently controlled incubation chambers, the platform enables true high throughput, small scale bioprocessing. Integrated mixing improves oxygen transfer and reduces gradient effects, while continuous DO and pH readouts provide real time culture insight across parallel, independently controlled environments.



1 Patented 96/24-well plate mixing

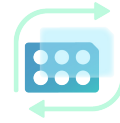


2 Real-time DO/pH monitoring




3 Four independent incubation chambers



S.NEST 2.0

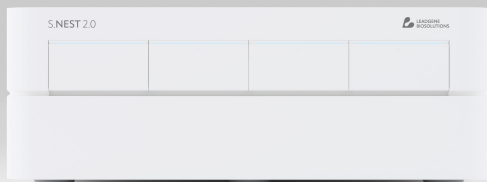

C.NEST 2.0

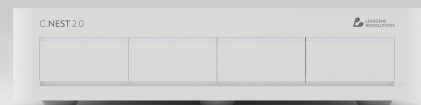

C.BIRD



# Patented Reciprocal Mixing Technology

S.NEST®, C.NEST®, and C.BIRD™ apply patented pneumatic mixing technology via a proprietary mixing lid, delivering precise, programmable mixing in standard commercial 96- and 24-well plates.


Gentle, adjustable mixing enables true beneficial agitation even at the 96-well format. Effective and reproducible mixing at the 24-well scale to achieve shaker-ready cell numbers.

Applied in CHO single-cell workflows, the platform enables faster, more consistent clone expansion and selection.

- ⑤ Uniform nutrient distribution
- ⑤ Higher oxygen transfer




Air-driven suction and expulsion enable efficient pneumatic mixing in 96- and 24-well plates, with adjustable mixing rates to match your culture conditions.

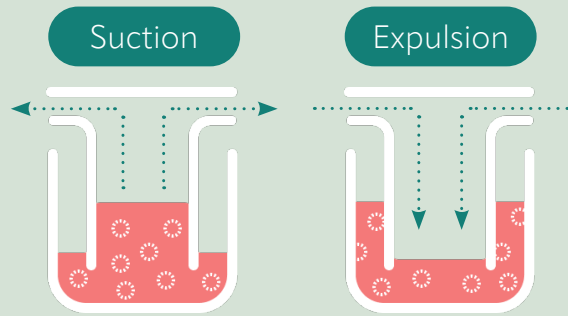


## Special feature - Suspension boost

Suspension Boost is a high energy mixing mode designed to support higher viable cell densities in 24-well plates. It is particularly useful for fed batch workflows and medium optimisation studies.

In CHO scale down models, this mode can increase achievable cell density by five to ten times compared to static conditions, depending on the application.

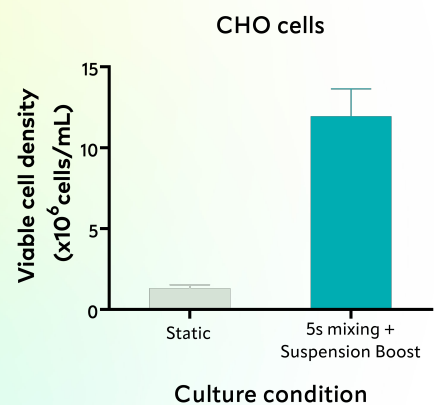

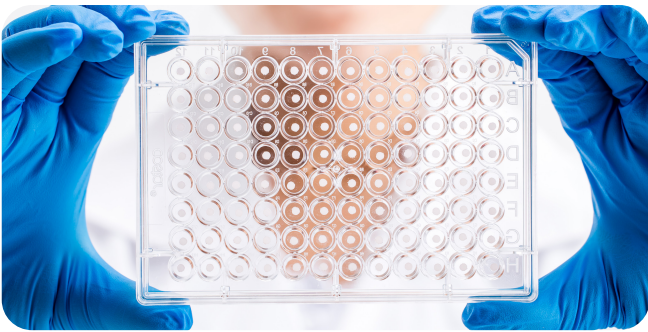


# Real-Time DO/pH Monitoring

Measurements are non-invasive, enabling continuous monitoring without disrupting cultures or increasing contamination risk.

↳ Monitor cell growth trends and metabolic activity in real time


↳ Collect measurements every 10 minutes across all wells

S.NEST measures DO and pH in every well using optical sensor spots integrated into standard commercial 96- and 24-well plates.

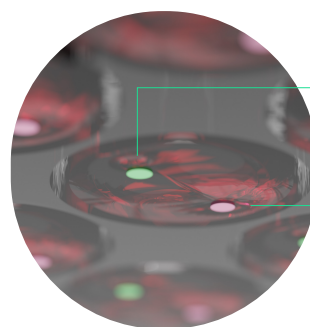


Dissolved oxygen  
sensor tag

pH level  
sensor tag

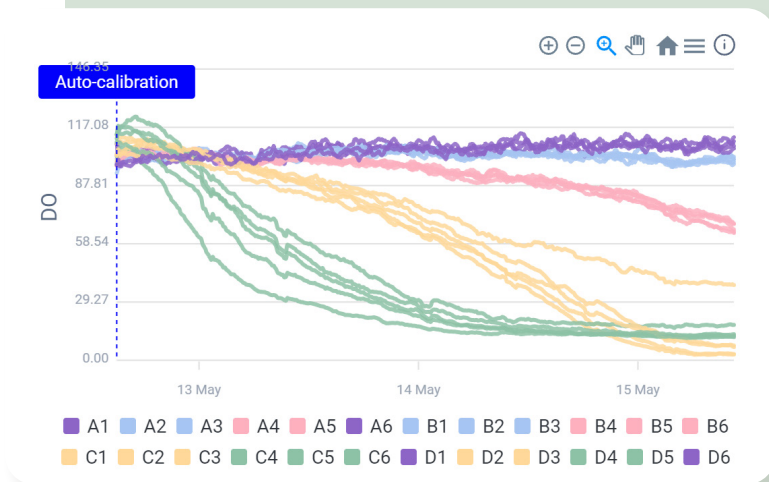

- Control
- $1 \times 10^5$  cells/mL
- $2 \times 10^5$  cells/mL
- $5 \times 10^5$  cells/mL
- $1 \times 10^6$  cells/mL

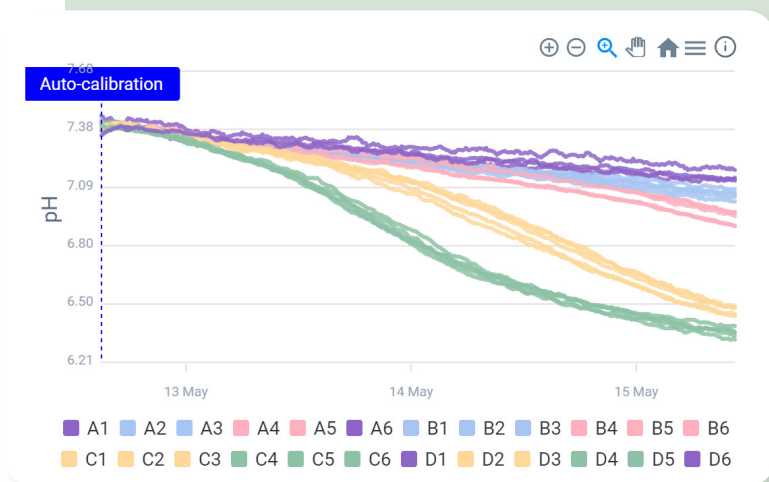

## Real-time indicators for clone selection and medium optimisation



### DO monitoring

Cell growth rate  
& Cell viability

↘ Cell growth profile and  
harvest time point



### pH monitoring

Lactate accumulation  
& Metabolic switch

↘ Feed and pH control  
time point


# Optimized In-Well Environment & Multi-Chamber Scalability

Cultures in the four chambers can be independently configured with different incubation settings, mixing patterns, and sensing options, enabling parallel experiments with greater flexibility.



## Four independently controlled incubation chambers

- Disposable water reservoir
- Temperature control
- CO<sub>2</sub> level control
- Humidity monitoring
- UV light sterilization

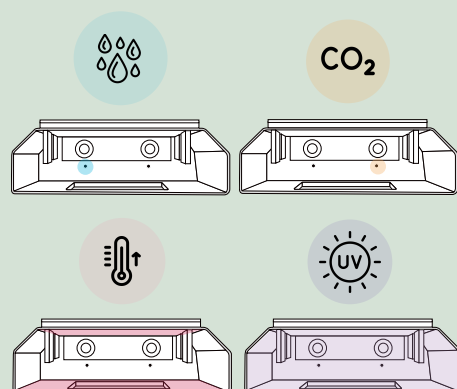



## Choose the system that fits your workflow

	<b>S.NEST 2.0</b>	<b>C.NEST 2.0</b>	<b>C.BIRD</b>
Mixing rate (Sec/cycle)	5–50 (±5%)	5–50 (±5%)	15-50 (±2%)
Suspension boost	Yes	Yes	N/A
Throughput	96 * 4 / 24 * 4	96 * 4 / 24 * 4	96 * 3 / 24 * 3
Medium Volume (µL/well)	1000 – 1600 150 – 200	1000 – 1600 150 – 200	1000 – 1600 150 – 200
Incubation chambers	4	4	N/A
Temperature Control (°C)	(RT+10) to 40	(RT+10) to 40	N/A
CO <sub>2</sub> Control (%)	1 – 20	1 – 20	N/A
Humidity Monitoring	Yes	Yes	N/A
UV light Sterilization	Yes	Yes	N/A
Data Output	DO / pH	N/A	N/A
Automation Compatibility	Yes	Yes	N/A
Workstation	Yes	Yes	N/A

# Accelerate CLD Workflows by 29–42%





## Accelerating Early Scale-Up in Cell Line Development

- Accelerates early scale-up by enabling faster cell expansion and a direct transition to larger formats.
- Reduces passaging by skipping intermediate 12- and 6-well steps.

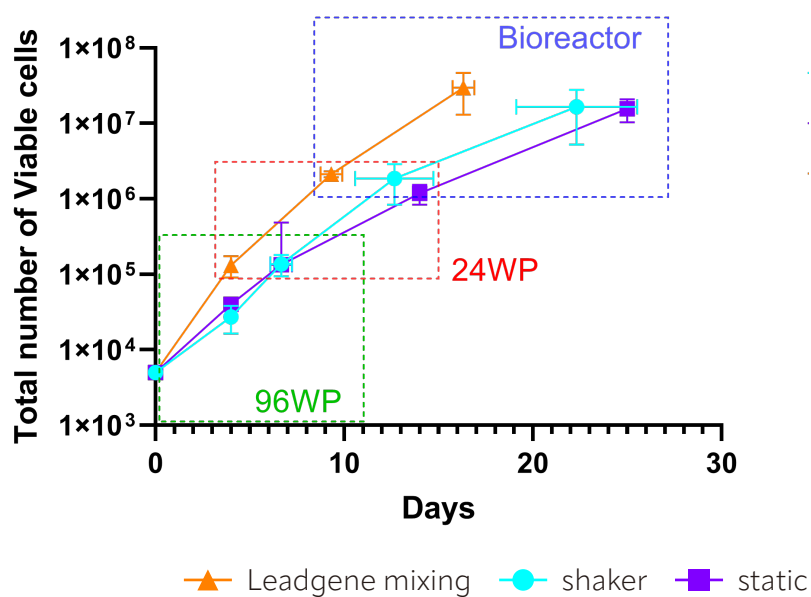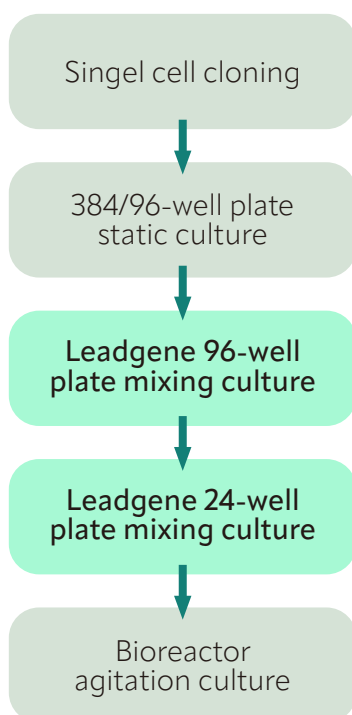

## Building Predictive Models for Process Development

- High-throughput scale-down models that improve clone ranking consistency with shake-flask results.
- Integrated DO/pH monitoring for real-time process insight.
- Supports data-driven and AI-assisted process optimization.

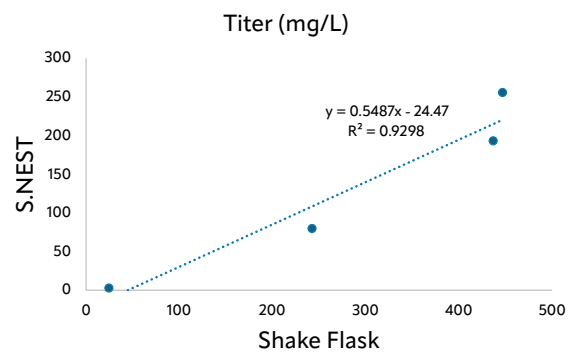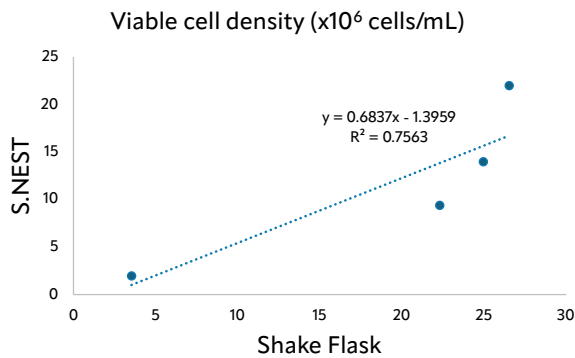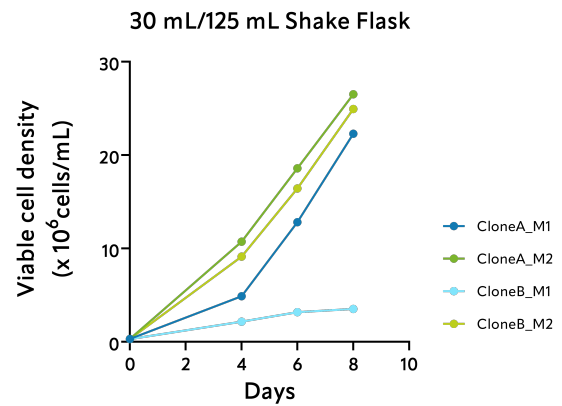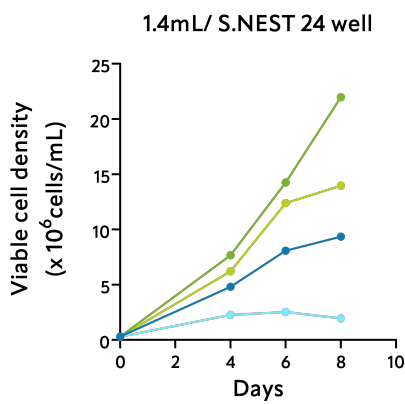

Accelerating Biology.  
Empowering Discovery.

LEADGENE BIOSOLUTIONS

[support@leadgene-bio.com](mailto:support@leadgene-bio.com) | [www.leadgene-biosolutions.com](http://www.leadgene-biosolutions.com)



©2025 Leadgene Biosolutions. All rights reserved. Duplication and/or reproduction of all or any portion of this document without the express written consent of Leadgene Biosolutions is strictly forbidden. Nothing contained herein shall constitute any warranty, express or implied, as to the performance of any products described herein. Any and all warranties applicable to any products are set forth in the applicable terms and conditions of sale accompanying the purchase of such product. Leadgene Biosolutions provides no warranty and hereby disclaims any and all warranties relating to the use of any third-party products or protocols described herein. The use of products described herein is subject to certain restrictions as set forth in the applicable terms and conditions of sale accompanying the purchase of such product. Leadgene Biosolutions may refer to the products or services offered by other companies by their brand name or company name solely for clarity and does not claim any rights to those third-party marks or names. Leadgene Biosolutions products may be covered by one or more patents. The use of products described herein is subject to Leadgene Biosolutions' terms and conditions of sale and such other terms that have been agreed to in writing between Leadgene Biosolutions and the user. All products and services described herein are intended FOR RESEARCH USE ONLY and NOT FOR USE IN DIAGNOSTIC PROCEDURES.

---